




## Mantelschürze BioClean-C

pure<sup>11</sup>-Nr.: 03207, Hersteller: BioClean

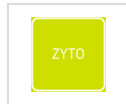

### Zusammenfassung

- Neue pure11-Artikelnummer (ab 01.07.2023): 1103207
- Material: Clean-Tough
- Zyto-geeignet
- Strapazierfähiges, leichtes Material
- Mantelschürze zum Binden
- Verstellbarer Halsabschluss
- Einzeln verpackt
- Geeignet bis zu ISO-Klasse 4

### Empfohlene Reinraumklassen

ISO	<input type="checkbox"/>	<input type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
GMP	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>

### Produktvarianten

**pure<sup>11</sup>-Nr.: 03207BS**

Farbe: Blau / Größe: S / Herst.-Nr.: BCAS / VE: 40 Stück

**pure<sup>11</sup>-Nr.: 03207BM**

Farbe: Blau / Größe: M / Herst.-Nr.: BCAS / VE: 40 Stück

**pure<sup>11</sup>-Nr.: 03207BL**

Farbe: Blau / Größe: L / Herst.-Nr.: BCAS / VE: 40 Stück

Quelle: <https://www.pure11.de/mantelschuerze-bioclean-c-0>

## About

The BioClean-C Chemotherapy Protective Apron with Sleeves provides protection against a range of cytotoxic drugs\*\*. Manufactured from lightweight low-linting CleanTough material, the apron features an adjustable neck fastening, tie tapes at the rear and ultrasonically bonded sleeve seams with protective tape.

## Specifications

**MATERIAL:**

BioClean-C CleanTough blue material

**COLOUR:**

Blue

**CONSTRUCTION:**

Ultrasonically bonded sleeve seams with protective tape & 100% polyester elasticated cuffs.

**CLASSIFICATION:**

PPE Cat 3 Type PB [6]

**CHARACTERISTICS:**

Low particulating

**COMPATIBILITY:**

ISO Class 4

## Features

- Tested against permeation standard ASTM F739-12
- Ultrasonically bonded seams with protective tape
- Tie tapes at rear
- Adjustable neck fastening
- Lightweight low-linting CleanTough material
- Silicone-free

EN 13034:2005  
+ A1:2009



TYPE PB[6]



**LIGHTWEIGHT & LOW-LINTING**

**TESTED AGAINST PERMEATION STANDARD ASTM F739-12**

**ULTRASONICALLY BONDED SEAMS**



Dr Isolator© represents BioClean's GMP compatible Chemo Safety Wear

## Quality Standards

- Manufactured in a facility holding ISO 9001:2015 and (EU)2016/425 Module D certifications
- Conforms to Category 3 - Personal Protective Equipment Regulation (EU) 2016/425
- Tested against standard EN ISO 13034:2005 + A1:2009 protective clothing against liquid chemicals (Type PB [6])
- Tested against standard EN 14325 protective clothing against chemicals
- Tested against permeation standard ASTM F739-12
- Processed in an NEBB certified ISO Class 4 environment
- Tested against standards: EN 530, Abrasion resistance; EN 6530, Resistance to penetration by liquids; EN 13934-1, Tensile strength; EN 9073-4, Trapezoidal tear resistance; EN 863, Puncture resistance; and EN 13935-2, Seam strength.

## Shelf Life & Storage

Five (5) years from date of manufacture. Store in a dry, cool place (<40°C) away from direct sunlight and fluorescent light.

## Performance Test Results – BioClean-C CleanTough Blue Material

TEST	RESULT		PERFORMANCE CLASS (EN 14325)
Abrasion Resistance	10 to 100 cycles		1
Trapezoidal Tear Resistance	MD 50.6 N		3
Trapezoidal Tear Resistance	CD 32.3 N		2
Tensile Strength	CD 47 N		1
Tensile Strength	MD 94 N		2
Puncture Resistance	10 N		1
Repellence to Liquids	30% H <sub>2</sub> SO <sub>4</sub>	99.0%	3
	10% NaOH	98.4%	3
	O-Xylene	91.4%	1
	Butan-1-ol	93.9%	3
Penetration by Liquids	30% H <sub>2</sub> SO <sub>4</sub>	0%	3
	10% NaOH	0%	3
	O-Xylene	0%	3
	Butan-1-ol	0%	3
Seam Strength <sup>1</sup>	74 N		2

<sup>1</sup>Seam not destroyed

## Particle Shedding Test Results

TEST	RESULT	CATEGORY (IEST-RP-CC003.3)
Particle Shedding (Helmke Drum Test)	<1700	Category I

## BioClean-C™ CleanTough Blue Material - ASTM F739-12 Test Method Results

Mean Breakthrough Time (MBT), Minutes

Breakthrough of the test chemical is deemed to have occurred when the permeation rate has reached 0.1 µg/cm<sup>2</sup>/min

CISPLATIN	CARMUSTINE	CYCLO-PHOSPHAMIDE	DOXORUBICIN HYDROCHLORIDE	5-FLUOROURACIL	METHOTREXATE	ETOPOSIDE	PACLITAXEL	THIOTEPA
>240	>240	>240	>240	>240	>240	<6	>240	55 (46,88,31)

Results achieved under controlled laboratory conditions, by accredited external testing laboratory.

### TO ORDER

Re-Order	Chest	Height	Size
BCAS-S	84-92cm (33"-36")	164-170cm (5'4"-5'6")	Small
BCAS-M	92-100cm (36"-39")	170-176cm (5'6"-5'9")	Medium
BCAS-L	100-108cm (39"-42")	176-182cm (5'9"-6'0")	Large
BCAS-XL	108-116cm (42"-45")	182-188cm (6'0"-6'2")	X-Large

### Packing

One piece per sealed inner PE bag; one inner bag per sealed outer PE bag; 40 outer bags per lined carton (40 pieces).

For additional information visit us at [www.bioclean.com](http://www.bioclean.com), or call us at +44 1638 663338

**Nitritex Ltd**  
Minton Enterprise Park  
Oaks Drive  
Newmarket, Suffolk  
CB8 7YY  
United Kingdom

BCAS/1906/PDS47

Neither this document nor any other statement made herein by or on behalf of Ansell should be construed as a warranty of merchantability or that any Ansell product is fit for a particular purpose. Ansell assumes no responsibility for the suitability or adequacy of an end user's selection of aprons/sleeve covers for a specific application.

\*\*Please see product validation pack or contact Ansell customer service for specific data on use of garments with cytotoxic drugs. Garments used for protection against such drugs must be selected specifically for the type of chemicals used.

Ansell, ® and TM are trademarks owned by Ansell Limited or one of its affiliates. © 2012 – 2019 Ansell Limited. All Rights Reserved.

