




Tupfer TOC

Zubehör-Set, 72 Flaschen und 72 Etiketten
pure¹¹-Nr.: 06381, Hersteller: ITW Texwipe

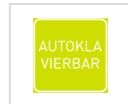

Zusammenfassung

- Neue pure11-Artikelnummer (ab 01.07.2023): 1106381
- Ersatzfläschchen und Etiketten
- Autoklavierbar
- Set besteht aus: 72 x 40 ml clear vials bonded septa caps (certified <10 µg/L TOC), 72 x blank vial labels

Empfohlene Reinraumklassen

ISO	<input type="checkbox"/>	<input type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
GMP	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>

Produktvarianten

pure¹¹-Nr.: 06381

Herst.-Nr.: TX3350 / VE: 1 Stück

Quelle: <https://www.pure11.de/tupfer-toc-2>

T E C H N I C A L D A T A S H E E T



TX3342 Bulk



TX714K Featured in **TX3340** TOC Cleaning Validation Kit



TX761K Featured in **TX3343** TOC Cleaning Validation Kit

Description

Texwipe's TOC Cleaning Validation Kits were designed to simplify sampling as part of a cleaning validation protocol. The swabs and vials are further cleaned using a proprietary process to make these kits excellent for use in Total Organic Carbon (TOC) analysis.

Lot coded for traceability and quality control.

Features & Benefits

Kit

- Organized, convenient and secure method for transporting swabs and vials from storage area to sampling sites to laboratory with minimal chance of contamination.
- Two size options available.

Swabs

- Certified Low TOC level (<50 µg/L, <50 ppb) ensures consistency in swab background contribution.
- Double layered, double knit polyester head entraps contaminants in knit structure during sampling. This allows maximum contaminant collection and release into the diluent to provide for excellent recovery rates.
- Notched, break-away handle allows the head to be placed into the vial with minimal handling and contamination.
- 100% polypropylene handle ensures no additional contaminants are introduced and provides excellent chemical resistance.

Vials

- Certified Low TOC level (<10 µg/L, <10 ppb) ensures consistency in vial background contribution.
- Prewashed vials have a one-piece cap and septa design for easy sampling and analysis.
- Polyethylene overcap protects septum surface from contamination until time of analysis.
- Sloped shoulder design minimizes headspace for reduced background TOC level.
- Endorsed for use by TOC instrumentation manufacturers.

Labels

- Cleanroom compatible for vial sample identification and traceability.

Container/Box

- Packed in a cleanroom compatible polypropylene box.

Applications

- Surface sampling for use in:
 - TOC analysis as part of a cleaning validation protocol.
 - As part of other cleaning validation protocols.
- Non-organic liquids sampling for contamination source investigations.


Products

See reverse for Physical and Contamination Characteristics.

Sterile Available

All Texwipe swabs available in sterile upon request.

Certificate of Analysis



Product Number: TX3342
 Lot Number: 133636
 Manufacture Date: 9/24/2013

CERTIFICATE OF ANALYSIS


Sorption			
Test	Specifications	Units	Result
Sorption Capacity	> 300	ul/swab	549

TOC			
Test	Specifications	Units	Result
Total Organic Carbon: Swabs	< 50	µg/L	46.7
Total Organic Carbon: Vial	< 10	µg/L	3.5

CERTIFICATE OF COMPLIANCE

ITW Texwipe certifies that the above product/material listed conforms to procedures and specifications according to ITW Texwipe Operating Procedures and Work Instructions.

ITW Texwipe certifies that the lot number provided under this Certificate has been inspected and tested, in accordance with ITW Texwipe specifications.

Approved By: 

Christopher R. Creasy

Date: **October 07, 2013**

Time: **12:53:44 PM EST**

QA-REC-7.5-002 REV.004 **Confidential**

ITW Texwipe
 1210 South Park Drive • Kernersville, NC 27284 • Tel: 800-839-9473

Instructions for Use

1. Please consult your written SOPs for sampling instructions.
2. Dip the swab in solution.
3. Sample the area.
4. Remove the vial cap.
5. Snap the swab head by placing the swab head into the vial with the notch touching the top of vial. Avoid touching the swab head.
6. Replace the cap.
7. Label the sample.

Products

Number	Description	Packaging
TX3340	Cleaning Validation Kit Components <ul style="list-style-type: none"> 12 40 mL clear vials bonded septa caps, certified <10 µg/L TOC 24 TX714K Low TOC Alpha® Sampling Swabs, certified <50µg/L TOC 12 blank vial labels 	18 kits/case
TX3342	Bulk Cleaning Validation Kit Components <ul style="list-style-type: none"> 72 40 mL clear vials bonded septa caps, certified <10 µg/L TOC 144 TX714K Low TOC Alpha® Sampling Swabs, certified <50µg/L TOC 72 blank vial labels 	1 kit/case
TX3343	Cleaning Validation Kit Components <ul style="list-style-type: none"> 12 40 mL clear vials bonded septa caps, certified <10 µg/L TOC 24 TX761K Low TOC Alpha® Swab with Long Handle, certified <50µg/L TOC 12 blank vial labels 	18 kits/case
TX3350	Low TOC Vials Kit Components <ul style="list-style-type: none"> 72 40 mL clear vials bonded septa caps, certified <10 µg/L TOC 72 blank vial labels 	1 kit/case

* Testing Method: TM2: Laboratory Testing for Swabs. Test method is available upon request. Values are typical, not representing specification limits.

Texwipe Regulatory Certificates

Texwipe is the only wiper company to be registered: ISO 9001 • ISO 13485 • ISO 14001 • OHSAS 18001